

Freeze yeast cells (make frozen stock)

Grow cells on YEP-D plates

Take cells by tooth pick, and put into 1ml of steriled 25%-glycerol solution.

(label clearly!)

Store at -80 degree.

or

Grow cells in YEP-D liquid media

Put 250ul of cell culture into screw cap tube

Add 250ul of steriled 50% glycerol solution.

Store at -80 degree.

Freeze e coli cells (make frozen stock)

Grow cells in LB-Amp liquid media

Put 250ul of cell culture into screw cap tube

Add 250ul of steriled 50% glycerol solution.

Store at -80 degree.